



Micropipet Protocol Card - Oxford

NEVER crank the micropipets above or below their ranges!!

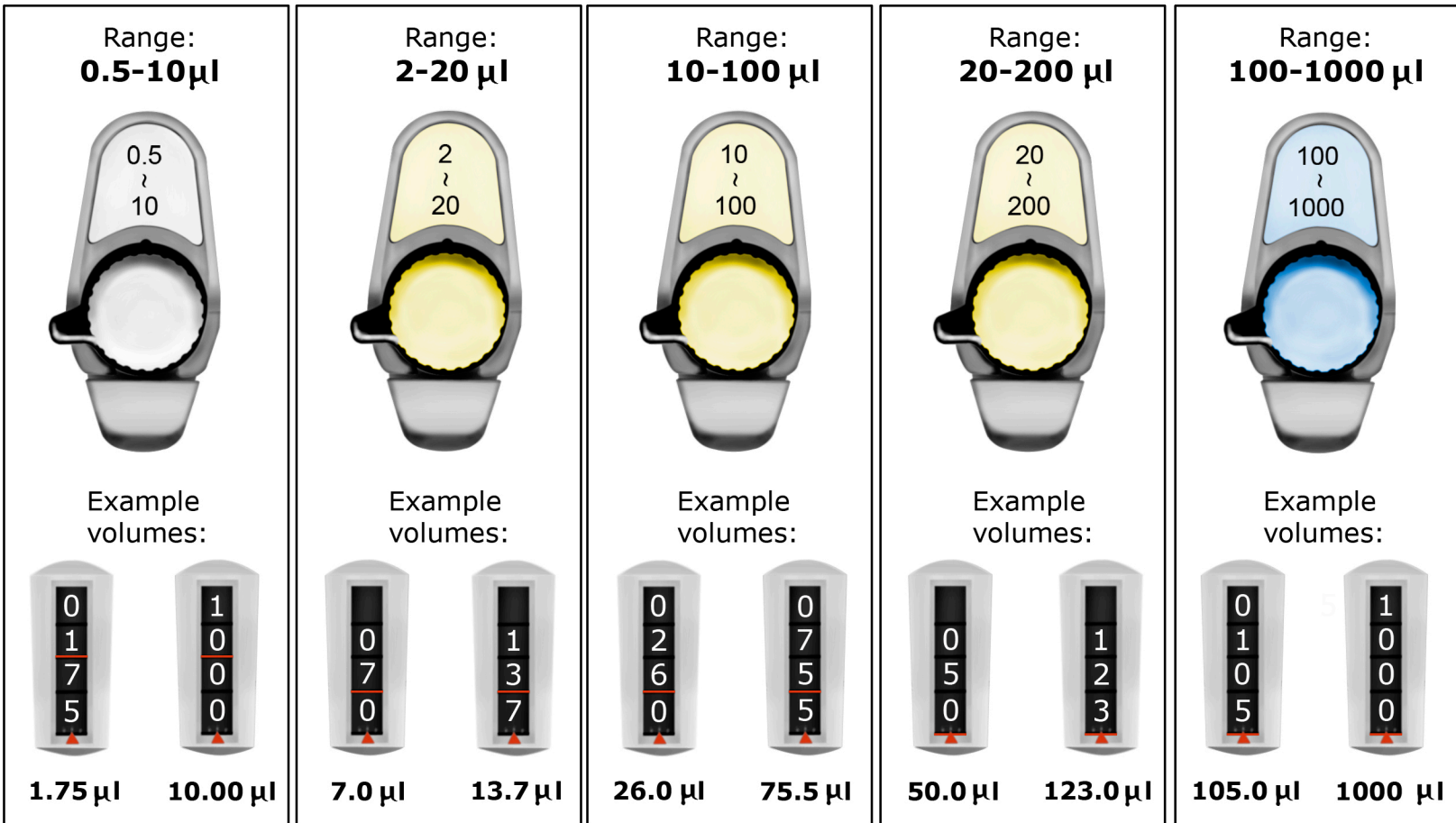


The red line indicates the decimal point.

Tip Selection Guide

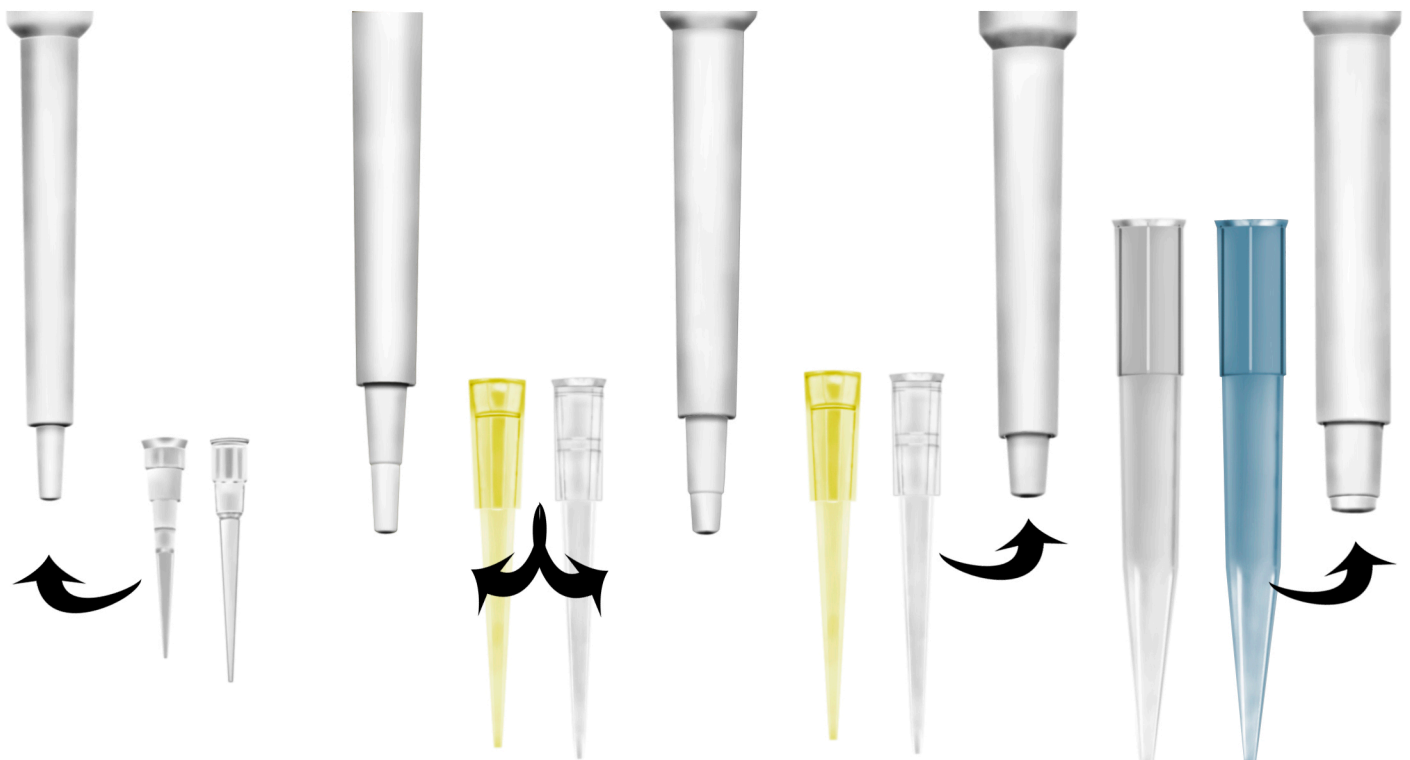

Always hold the micropipet in a **vertical position** when there is fluid in the tip.

Micropipet Protocol Card - Oxford

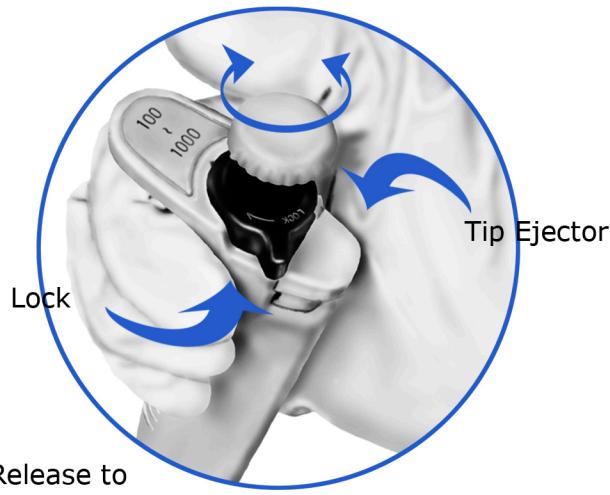

- To set the volume:
- 1) Unlock the knob
 - 2) Turn the knob to increase/decrease amount drawn up
 - 3) Lock knob

Push down to first stop

Release to draw in sample

Move to new microtube

Push down to first stop

Continue down to second stop

Eject Tip

