



A. Create a New User

1. Push "User" (F5)
2. Push "New" (F2)
3. Enter name by toggling with the arrow keys. Press enter to enter each letter or symbol selected as a name.
4. When the name is complete, press "Accept" (F1)

B. Create a Program

1. Press "User" (F5)
2. Press "Accept" (F1) after highlighting the correct user with the arrow toggle
3. Press "Create" (F2)
4. Edit the existing template program by using the arrow toggle and entering different values where desired. You do not need to hit the enter key. Move to the next value when you are done entering the new value.
5. When finished with the program, push "Store"
6. Push "Method" (F3) and name the method. Make sure to clear "exp000" or it will be incorporated into your name.
7. Push "Accept"(F1) when you have finished naming the program.
8. The machine will return to the main menu.

C. Run a Program

1. Push "User" (F5)
2. When the correct user name is highlighted, push "Accept" (F1)
3. Press "Run" (F1)
4. If there is more than one program stored under the user name, use the arrow toggle to highlight the desired program, then press "Start" (F1).
5. Enter the reaction volume and leave the ramp speed at 9600.
6. Press "Start" (F1) and the program will begin.
7. You may also view the desired program after you highlight it, but before you press start, by pressing "View" (F2)